




Here is the message sent by windows when trying to open the IOlab.exe application.

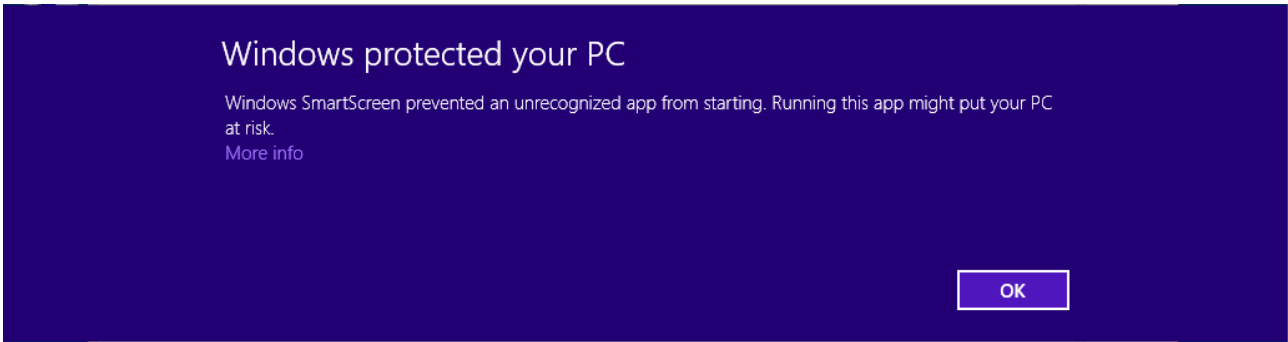


Clicking on more info, it is possible run anyway the app

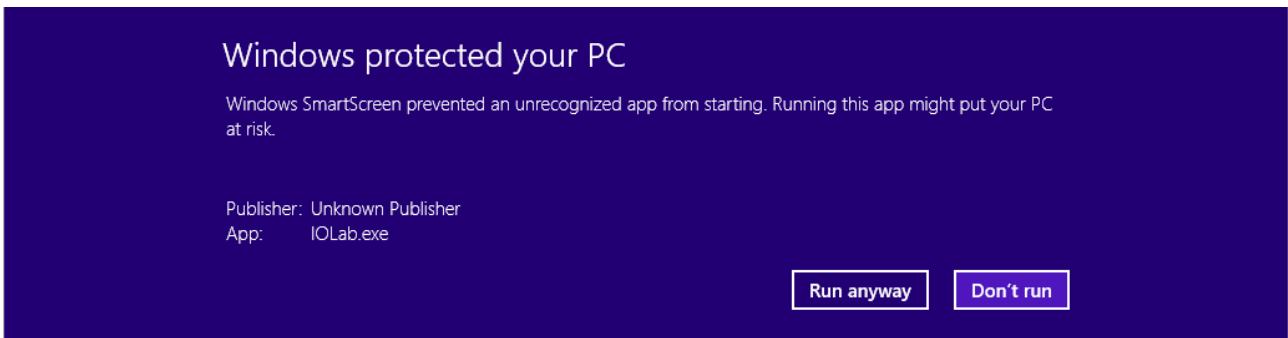

Then, windows complains it cannot open it because the file disappeared (Norton put it in quarantine)

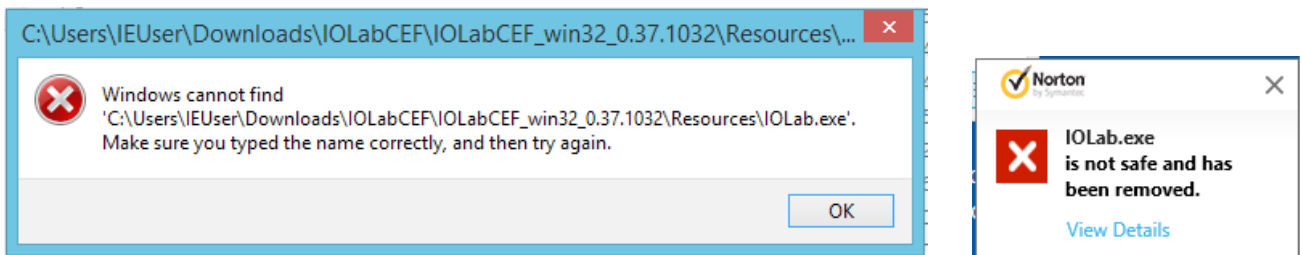

Open Norton and go to the History section

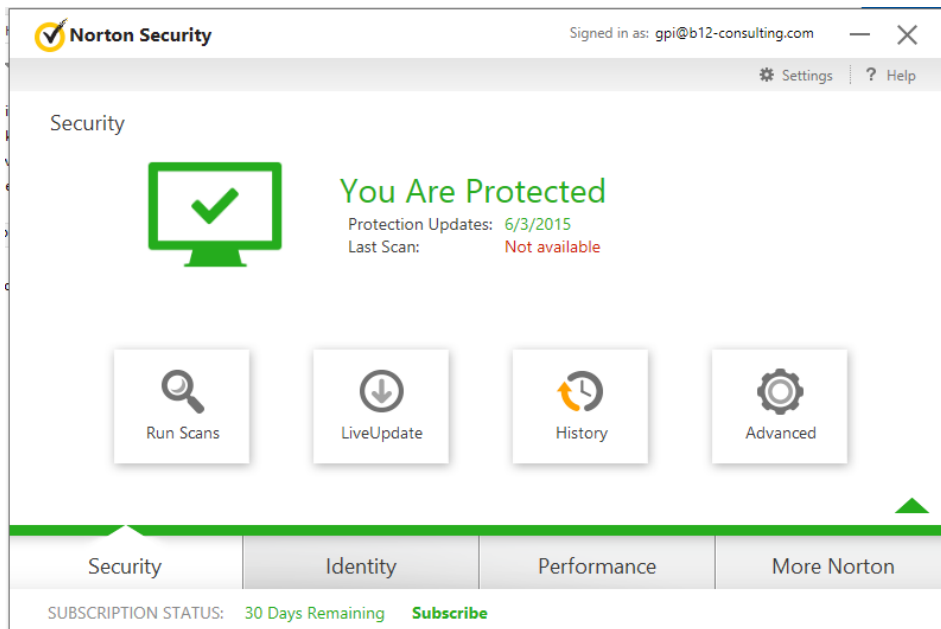

Security History ?

Show Recent History ↻ Quick Search Go

Severity	Activity	Status	Date & Time
● Medium	Download Insight analyzed IOlab.exe	Removed	9/2/2015 4:23:59 PM
● Medium	iolab.exe (WS.Reputation.1) detected by Download Insight	Quarantined	9/2/2015 4:23:59 PM
● Info	Firewall rules updated	Detected	9/2/2015 4:22:51 PM
● Info	You allowed Windows host process (Rundll32) to access your network resources.	Allowed	9/2/2015 4:22:48 PM
● Info	IPS Detection Statistical	Submitted	9/2/2015

Go to Page: Go Page 1 of 1

Details

Recommended Action
Resolved - No Action Required
[Restore](#) [Options](#)

! iolab.exe contained threat **WS.Reputation.1**

Risk
Medium

Origin
Not Available

Activity
Threat Actions performed: 1

[More Options](#)

[Import](#) [Export](#) Clear List Close

Then, select the entry for iolab with status Quarantined.
On the right panel, it is possible to restore the application...